


Microscope Stand Boom-Stand


HOW TO ASSEMBLE BOOM STAND – MDHA485S TUTORIAL

- Hold the black sleeve in place with the thermal head and tighten the side handle.

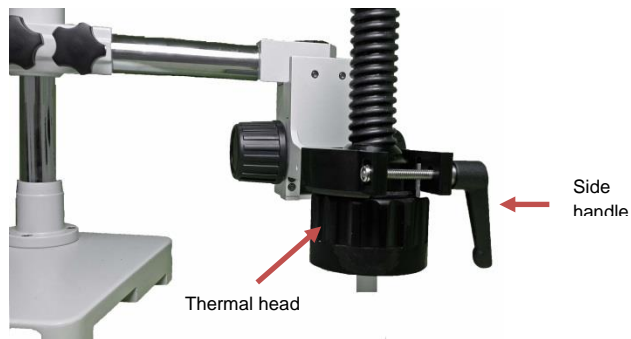

Fine tuning Z-Axis

- Rotate the Z axis **hand wheel's tension** to fine tune the plunger contact with the device.

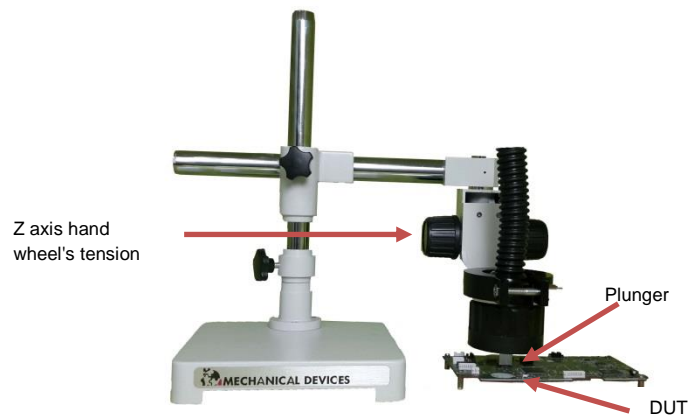

www.mechanical-devices.com

Package layout

- The Boom Stand is packed with parts laid out as shown in figure 1.


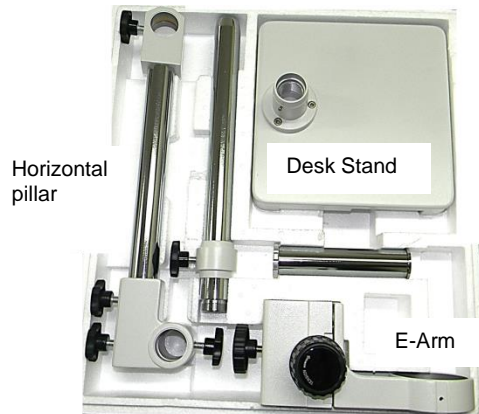



Fig. 1

Assembly

- Place the stand on top of the table and attached the bar tightly to place, with the assistance of small screw for stabilizing.

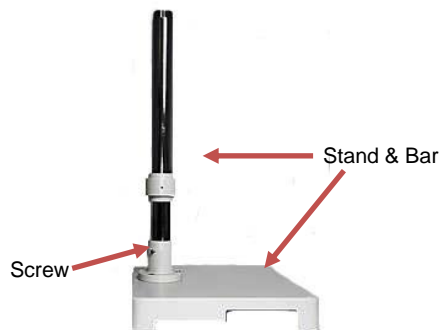

- Slide the **horizontal pillar** with cross block into the vertical post.

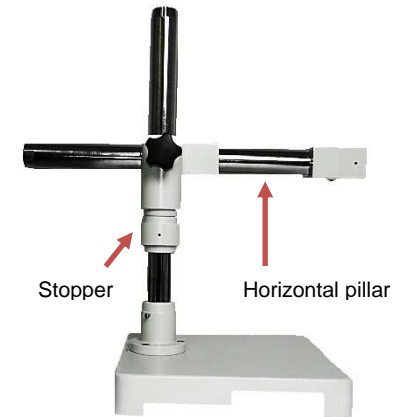

- Insert the e-arm into the end of the horizontal bar and tighten using lock wheel
- E-arm can be inserted from the TOP, or BOTTOM of the horizontal bar. Set e-arm up at various angles and tighten the lock lever for perfect oblique viewing.

